

Assembly Instructions

Part Number: EL-1

Product: EzyLok Kit


Standard accessory for all glass nebulizers.

1. P/N 70-808-0735/2: 1 x EzyLok 6mm Hose Adaptor
2. P/N QSM-4: EzyLok Connector for 4mm Tubing
3. P/N SNP-1/6: 1 x Hose Clip 6mm


Note: To release or connect the EzyLok Connector, press down firmly on the blue button.



Using the EzyLok Kit with EzyLok Arm Nebulizers

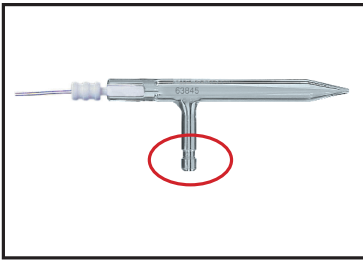


The EzyLok Arm

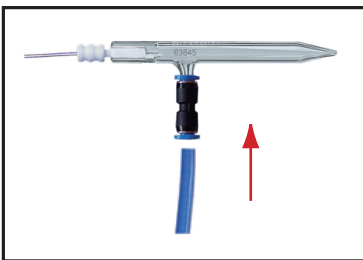


Step 2 (Option A):

If using Standard 4mm Tubing. Insert the tubing into the EzyLok Connector.

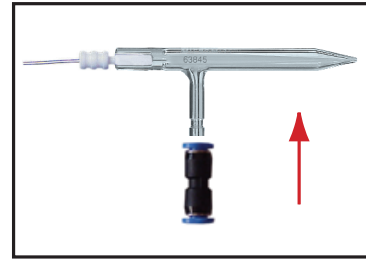

Step 1:

Connect the EzyLok Connector to the Nebulizer.

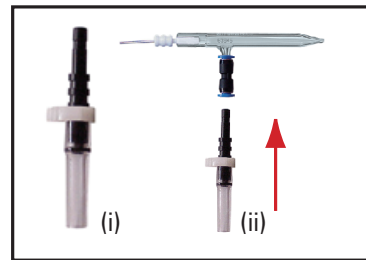

Step 2 (Option B):

If using other tubing. Secure tubing to Hose Adaptor with Hose Clip (i). The Clip should face the back of the nebulizer. Insert Hose Adaptor into Connector (ii).

Important Note: Before applying gas pressure, apply a slight amount of pulling pressure between the Nebulizer and Hose to ensure they are securely locked.



GLASS EXPANSION

Quality By Design

International

15 Batman Street

West Melbourne VIC 3003

Australia

A.B.N. 35 006 560 032

Telephone: +61 3 9320 1111

Toll Free (Aust): 1800 777 638

Facsimile: +61 3 9320 1112

Email: enquiries@geicp.com

www.geicp.com

Americas

4 Barlows Landing Road

Unit #2

Pocasset, MA 02559

USA

Telephone: 508 563 1800

Toll Free: 800 208 0097

Facsimile: 508 563 1802

Email: geusa@geicp.com